



# Tissue I.D.™

Catalog #6730

---

---

## Intended Use

Tissue I.D.™ is a dye designed to stain tissues for visualization after processing and to facilitate orientation at the embedding station.

---

---

## General Information

Tissue I.D.™ is used to stain tissues to help orient the specimen for embedding. Tissue I.D.™ is used at the grossing station and can be applied to fresh or fixed tissues. Tissue I.D.™ survives processing and lightly stains small biopsy specimens. The stained surface of the small specimens facilitates orientation of the tissues for embedding. The mucosal surface of the biopsy specimen is more easily identified with Tissue I.D.™

---

---

## Packaging

Catalog #	Volume
#6730	60mL Dropper Bottle
#6731	120mL Dropper Bottle x 4/cs

---

---

## Staining Procedure

Fresh or fixed tissue may be stained with Tissue I.D.™

1. With an applicator, dropper, or brush, apply Tissue I.D.™ to specimen.
2. Place the specimen in the desired fixative prior to placing on the processor.
3. The Tissue I.D. will survive processing and leave a lightly stained tissue surface to facilitate proper orientation for embedding.